



# SANIHOL<sup>®</sup> 70

## 70% Denatured Ethanol Solution

### Technical Data Sheet

#### - BRIEF DESCRIPTION -

*SANIhol is a 70% (v/v) denatured ethanol solution that is filtered to 0.2 microns. SANIhol can be used to clean surfaces in microbiology labs and tissue culture labs. It is also effective in cleaning surfaces and process equipment in pharmaceutical, biotechnology, and medical device, cosmetic and production areas. The alcohol solution evaporates away completely so no rinse is required.*

#### - USE PROTOCOL -

Use in well-ventilated area. Spray or wipe solution on surface to be cleaned until it is completely covered. Allow to air dry. When using trigger spray, point away from eyes and skin. Hold bottle six to eight inches from surface to be treated.

#### - INGREDIENTS-

63.4% Ethyl Alcohol  
3.2% Methyl Alcohol  
3.4% Isopropyl Alcohol  
30% USP Purified Water % by volume

#### - SHELF LIFE AND STORAGE -

The shelf life is 2 years from the date of manufacture. The expiration date can be found on the product. Do not store at temperatures above 120°F.

#### - PRECAUTIONARY STATEMENT -

**Flammable** - do not come in contact with flame or extreme heat. Avoid eye and skin contact and ingestion.

#### - PACKAGING -

**16 oz. Pump Spray / 12 per case**  
**1 Gallon / 4 per case**

#### - CATALOG NUMBERS -

<u>Manufacturer</u>	<u>Item</u>	<u>Cat. No.</u>
Decon Labs, Inc.	16 oz.	8616
	1 Gallon	8601

**For Additional Technical Information call  
Decon Customer Service  
(800) 332-6647, or  
See our website at [www.deconlabs.com](http://www.deconlabs.com)**

®Decon Labs, Inc.

Revised 9/13/2013