

KIMTECH PURE* A5 Sterile Cleanroom Apparel with CLEAN-DON* Technology

Gowning Procedure

BEFORE GOWNING

Step 1: (Pre-Entry) Don Hair Net and Shoe Covers after removing all jewelry and cosmetics.

Step 2: (Gowning) Wash hands and gown first pair of sterile gloves. Sanitize gloves after gowning each article if required.

Step 3: Apply mask and hood assuring a snug fit.

Step 4: Open vacuum-packed apparel.
Tear at notched edge.



Step 5: Grasp the blue line.
Located on the inside middle back.



Step 6: Gently unfold coverall utilizing blue indicator line. Arms and legs are pre-drawn and snapped in place. Garment is already folded inside-out and unzipped.

BEGIN GOWNING

Snap allow gathered-up arms and legs to expand during gowning



Step 7: Hold garment at waist



Step 8: Put one leg in and point toe through opening until snap releases



Step 9: Do the same with the other leg



Step 10: Insert one arm and extend until snap releases



Step 11: Do the same with the other arm



Step 12: Slip thumbs through thumb loops



Final Step: Cross legs and zip up coverall



After A5 Sterile Cleanroom Apparel is Gowned:

Add boot covers



Complete gowning by adding goggles and a second pair of sterile gloves.