

Kimtech Pure* A5 Sterile Cleanroom Apparel Gowning Procedure

Before Gowning

Step 1

(Pre-Entry) Don Hair Net and Shoe Covers after removing all jewelry and cosmetics.

Step 2

(Gowning) Wash hands and gown first pair of sterile gloves. Sanitize gloves after gowning each article if required.

Step 3

Apply mask and hood assuring a snug fit.

Step 4

Open vacuum-packed apparel. Tear at notched edge.

Step 5

Grasp the blue line. Located on the inside middle back.



Step 6

Gently unfold coverall utilizing blue indicator line. Arms and legs are pre-drawn and snapped in place. Garment is already folded inside-out and unzipped.



Snap fasteners allow gathered-up arms and legs to expand during gowning.

Begin Gowning

Step 7

Hold garment at waist.



Step 8

Put one leg in and point toe through opening until snap releases.



Step 9

Do the same with the other leg.



Step 10

Insert one arm and extend until snap releases.



Step 11

Do the same with the other arm.



Step 12

Slip thumbs through thumb loops.



Final Step

Cross legs and zip up coverall.



After A5 Sterile Cleanroom Apparel is gowned

- Add boot covers.
- Complete gowning by adding goggles and a second pair of sterile gloves.



Exceptional
Workplaces®

For more information, contact your local Kimberly-Clark Professional® Sales Representative at 1-800-255-6401 or visit www.kimtech.com.

KIMTECH
PURE* BRAND