

# KIMTECH PURE\* M3 Sterile Pouch Face Mask Donning Procedure

The Science of Comfortable Protection.



## Pre Hood Donning Procedure

### Step 1

Carefully open the protective packaging.



### Step 2

While holding the mask through the protective packaging, reach in and separate the head band straps.



### Step 3

Pull bands over the head and secure while holding the gap guard to keep mask in place.



### Step 4

Pinch nose wire into place.



### Step 5

Inhale and exhale to check for secure fit.



### Step 6

Put hood on over the mask

